
## 10720 : STAGE-TOP INCUBATOR FOR MÄRZHÄUSER STAGES WITH 160x110MM INSERTS

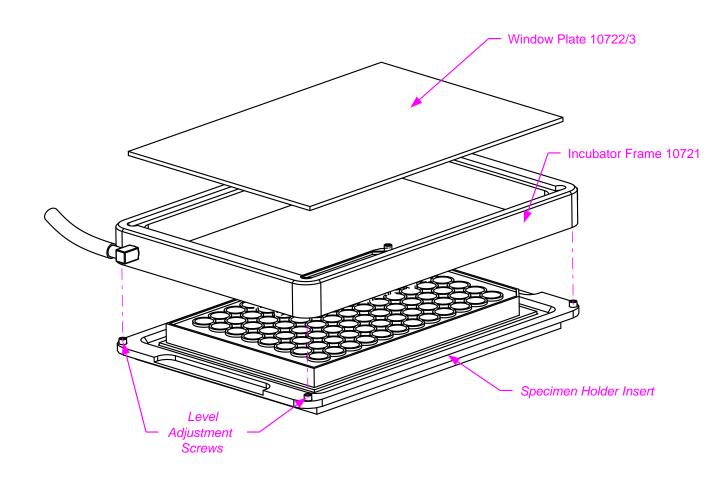

Insert a window plate into the the incubator frame 10721. Use the glass window plate 10723 if you plan to use contrast modes that require a polarization-maintaining material (e.g. interference contrast). Use the unbreakable polycarbonate window plate 10722 otherwise.

Adjust the four level adjustment set screws of the specimen holder insert to protrude from the surface. (Longer set screws can be used if they are needed for their original purpose also.

Connect the outlet of the humidifer column to one of the ports on the sides of the incubator frame as shown. Use the port on the opposite side if re-criculation of the gas is desired.

Place the incubator on the specimen holder insert as shown.