

13197 : STAGE-TOP INCUBATOR SEMI-OPEN FOR LEICA SUPER-Z GALVO STAGE WITH ROTATING INSERT AND S28 CONDENSER

Preparations:

- » Place a (dummy) dish or chamber container on the rotating insert of the Leica Super-Z galvo stage and close any gaps remaining in the insert around the dish as far as possible.
- » Connect the inlets of the incubator frame to the outlet of the humidifier column using flexible tubing and the Y-connector supplied.

Operation:

- » Move the galvo stage to the lowest position.
- » Place the incubator frame on the rotating insert of the Leica Super-Z galvo stage.
- » Tilt the lamp pillar, slide the incubator cover assembly over the S28 condenser, and put the pillar back in the vertical position.
- » Adjust the height of the incubator cover assembly so that it just touches the incubator frame.
- » Adjust the position of the cover disc and the cover ring to allow access to the sample when needed.

