
## 15035: STAGE-TOP INCUBATOR CLOSED 160x110x18mm

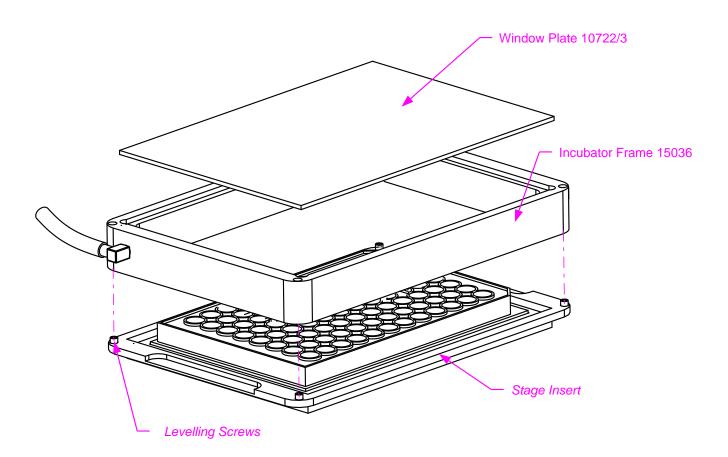

Insert a window plate into the the incubator frame 15036. Use the glass window plate 10723 if you plan to image in transillumination modes (e.g. interference contrast). Use the unbreakable polycarbonate window plate 10722 otherwise.

Adjust the four levelling screws of the stage insert so they protrude slightly from the top surface. (Longer set screws can be used if the levelling screws are needed for their original purpose also.

Connect the outlet of the humidifer column to one of the ports on the sides of the incubator frame as shown. Use the port on the opposite side if re-criculation of the gas is desired.

Place the incubator on the specimen holder insert as shown.